

Method Number: TM 004

Updated: 16/03/2022

Issue Number: 25

Page 1 of 1

**Method Summary****Solvent (Dichloromethane or Cyclohexane) Extractable Material from Soils****Scope and Range**

This method details the solvent (dichloromethane or cyclohexane) extraction of soil samples by Soxtherm to quantify the extractable organic material present.

Calibration Range: 100 mg/kg to 10⁶mg/kg

Sample Range: LOD - 10⁶mg/kg

References

none

Principle

A known mass of dried and crushed soil sample is added to a Soxtherm thimble. The soil is heated in solvent (DCM or cyclohexane) under reflux. The solvent is then evaporated leaving a residue. The residue is weighed and related back to the initial mass of sample to give the concentration of extractable matter in the sample.

Interferences

Solvents, reagents glassware and other sample processing hardware may yield artefacts and/or interferences under the conditions of analysis. This is undertaken by analysis of extracted blanks.